

# Delvotest® is easy to use



- Switch on incubator and leave to reach 64°C
  - 1 Use representative sample of milk
  - 2 Remove ampoule from the package – be careful not to damage the foil of the remaining ampoules (Always store spare ampoules in the fridge between 4°C and 8°C)
  - 3 Open and identify the ampoule
  - 4 Collect milk into the pipette
  - 5 Pipette the milk into the ampoule
  - 6 Verify the temperature (64°C +/- 2°C) of the incubator before placing the ampoule in
  - 7 Place the ampoule into the incubator immediately if it is on 64°C and press start to run the control time.
- Read the colour\* of the solid agar in the ampoules after the required incubation period\*\*



**Milk ok**

**Milk not ok**

\* The colours are indicative

\*\* Before using Delvotest® make sure to read the instructions for use

\*not listed, legislation regulation normally restricts the use: not to be used for lactating animals.

Antibiotic class	Antibiotic	Examples of branded products containing the antibiotic ingredient alone or in combination.	MRL (EU)	Delvotest® SP-NT (Unit: ppb)	Delvotest® T (Unit: ppb)
				Ampoules	Ampoules
Penicillins	Amoxicillin	Duphamox LA; Synulox	4	2	2
	Ampicillin	Duphacillin	4	2	2
	Penicillin G	Duphopen; TetraDelta; Ubro Red DC	4	2	3
	Cloxacillin	Kloxerate, Orbanin	30	12	10
	Oxacillin		30	4	3
	Nafcillin		30		3
Tetracyclines	Oxytetracycline	Engemycin LA; Terrramycin LA	100	300	100
	Chlortetracycline		100	300	150
	Tetracycline		100	300	80
	Doxycycline		*	150	50
Sulfonamides	Sulfamethazine	Duphatrim IS; Norodine 24; Lactatrim MC	100	75	125
	Sulfathiazole		100	40	30
	Sulfadimethoxine		100	50	40
	Sulfadiazine	Diatrim	100	65	55
	Sulfadoxine		100		80
Macrolides	Tilmicosin	Mycotil	50	30	60
	Tylosin A	Tylan 200	50	35	35
	Erythromycin A		40	90	160
	Lincomycin		150	170	
	Rifaximin		60	50	
	Spiramycin		200		1500
Aminoglycosides	Neomycin B	Multiject IMM; TetraDelta	1500	115	140
	Gentamycin		100	90	100
	Kanamycin	Ubrolexin	150	1700	
	Streptomycin		200		700
	DHStreptomycin	TetraDelta	200	700	700
	Spectinomycin		200	2000	
Cephalosporins	Cephapirin	Mastiplan; Metricure	60	2	5
	Desacetylcephapirin		60		2
	Ceftiofur	Exenel (if used as directed, nil milk withdrawal)	100	20	20
	Desfuroylceftiofur		100		45
	Cefalonium	Cepritect	20		5
	Cefoperazone	Pathocef	50	30	20
	Cefalexin	Ubrolexin	100	45	30
	Cefquinome	Cephaguard; Cobactan MC	20	65	50
	Cefazolin		50		3
Others	Lincomycin		150		275
	Chloramphenicol		*	3000	4000
	Trimethoprim	Norodine, Lactatrim MC, Diatrim	50	160	110
	Rifaximin		60		40
	Pirlimycin		100		300
	Dapsone		*	2	10
	Clavulanic Acid		200		700

